

Replaceable Stopcock Assembly Instructions

GP1089, GP1090, and GP1083



Introduction

This procedure is for the assembly of the buret body and replaceable parts for the following Flinn products; GP1089, 25-mL buret, GP1090, 50-mL buret, and GP1083, 100-mL buret.

Procedure

1. Remove the split white nut from the stopcock assembly.

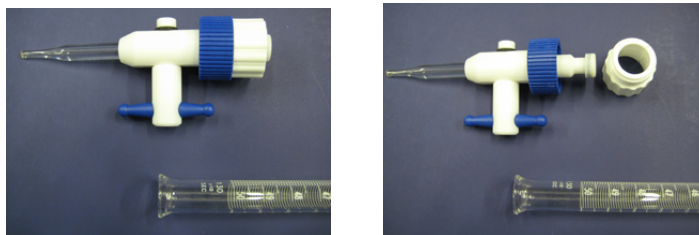


Figure 1.

2. Slip the split white nut onto the buret body making sure the threads face the bottom of the buret body (see Figure 2).



Figure 2.

3. Fully insert the stopcock assembly into the bottom of the buret body.

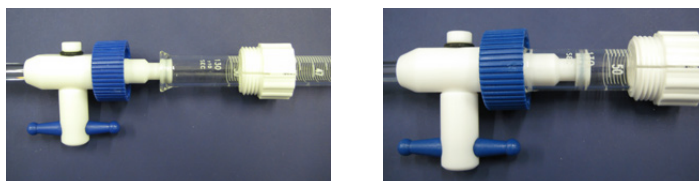


Figure 3.

4. Thread the split white nut into the blue nut of the stopcock assembly. Continue threading the nut until the fit is snug.



Figure 4.